

Semba Octave™ Chromatography System



User Manual Version 7.0

Rev 072809



Semba Biosciences, Inc.
5520 Nobel Drive, Suite 150, Madison, WI 53711
Phone: 608-310-4457 Fax: 608-310-4456




www.sembabio.com

CONTENTS

1.	ABOUT THE SYSTEM	5
1.1.	Description	5
1.2.	Supplied Components.....	5
1.3.	Components Not Supplied.....	7
1.3.1	Required Components.....	7
1.3.2	Optional Components.....	7
1.4.	Overall Specifications	7
1.5.	Pump Module	7
1.5.1	Pump Features	7
1.5.2	Pump Specifications	8
2.	INSTALLATION	9
2.1.	Precautions	9
2.2.	Location and Setup.....	9
2.3.	Unpacking and Inspection	9
2.4.	Location/Environment.....	9
2.5.	Pneumatic Connection	10
2.6.	Electrical Connections	10
2.7.	Inlet and Outlet Connections.....	12
2.7.1	Solvent Supply to Pumps.....	12
2.7.2	Pump Module Connections	12
2.7.3	Internal Inlet and Outlet Lines.....	12
2.7.4	Column Connections	13
2.7.5	Fittings	13
2.7.6	Pump Designations and Configurations.....	13
2.8.	Column “Jumper” Connections	14
3.	PUMP PREPARATION	16
3.1.	Solvent Preparation.....	16
3.1.1	Solvent Out-gassing and Sparging.....	16
3.1.2	Cavitation.....	16
3.1.3	Filtration	16
3.2.	Mobile Phase Reservoirs	17
3.3.	Self-Flush Solution	17
3.4.	Self-Flush System Set-up.....	17
3.5.	Priming the Pump Heads	17
3.6.	Long Term Pressure Calibration Accuracy.....	18
3.7.	Preparation of Pumps for Storage	18
4.	POWERING UP AND SHUTTING DOWN THE SYSTEM	19
4.1.	Powering Up the System.....	19
4.2.	Shutting Down the System	20

5.	SEMBAPRO™ SOFTWARE	21
5.1.	Starting the Program	21
5.1.1	Installation.....	21
5.2.	Using the Program	21
5.2.1	Opening Screen: the Current Valve State Window	21
5.2.2	Valve Names and Operation	22
5.2.3	Pump Control Window	23
5.3.	Using Scripts	24
5.3.1	Script Window	24
5.3.2	Creating a Script	25
5.3.3	Running a Script	26
6.	CONNECTING COLUMNS	28
6.1.	Purging the System	28
6.2.	Column and Back Pressure Considerations	28
6.3.	Connecting Columns	29
6.3.1	Column Rack	29
6.3.2	Connection to Valve Block Ports using ShapeShifter™ Adapters	29
6.3.3	Column Connection Procedure	29
7.	PERFORMING CHROMATOGRAPHY	33
7.1.	Running a Single Column	33
7.2.	Running Multiple Columns in Tandem	33
7.3.	Optimizing Chromatography Methods	33
7.4.	Classical SMBC Process	35
7.4.1	Designing SMBC Methods with the Octave System	36
7.5.	Step SMBC	37
7.5.1	Step Chromatography Configuration	37
7.5.2	Determining parameters for protein affinity purification in Step SMBC	38
7.5.3	Example Step SMBC method	39
7.5.4	Other Applications for Step SMBC	40
7.6.	Isocratic SMBC	40
7.6.1	“Isocratic Affinity” SMBC Purification of Proteins	41
7.6.2	Practical Considerations for Isocratic SMB-IMAC	42
7.6.3	Mathematical Model for Isocratic SMBC	42
7.6.4	Example: Chiral Separation Method Development	44
7.6.5	Optimizing SMBC Conditions	46
8.	SMBC PROTOCOLS FOR PROTEIN PURIFICATION	48
8.1.	Step SMBC for affinity purification	48
8.2.	Isocratic SMBC for affinity purification	50
9.	REFERENCES	52
10.	APPENDIX 1: MANUAL PUMP OPERATION	53
10.1.	Front Panel Controls and Indicators	53
10.1.1	Digital Display.....	53
10.1.2	Keypad.....	53

10.1.3	Status LEDs.....	53
10.1.4	Power-up Configuration.....	54
10.2.	Theory of Operation.....	54
10.2.1	Pump Cycle	54
10.2.2	Electronic Control	54
11.	APPENDIX 2: PUMP MAINTENANCE	56
11.1.	Mobile Phase Solutions.....	56
11.2.	Changing Pump Heads.....	56
11.2.1	Removing the Pump Head	56
11.2.2	Cleaning the Pump Head Assembly.....	57
11.2.3	Replacing Piston Seals.....	58
11.2.4	Changing the Piston	59
11.2.5	Replacing the Pump Head	59
11.3.	Conditioning New Seals.....	59
11.4.	Check Valve Cleaning.....	60
11.5.	Cleaning the Pump	60
11.6.	Cleaning the Cabinet	60
11.7.	Lubrication	60
11.8.	Fuse Replacement	60
11.9.	Battery Replacement	61
11.10.	Pump Troubleshooting	62
12.	APPENDIX 3: PREPARATION OF PROTEIN SAMPLES FROM BACTERIAL CELL CULTURES.....	63
12.1.	Cell culture and induction of target protein expression (pET vectors)	63
12.1.1	Autoinduction shake-flask protocol.....	63
12.1.2	IPTG induction shake-flask protocol.....	63
12.2.	Bacterial cell lysis and preparation of the soluble fraction (Feed).....	64
12.2.1	Cell lysis by detergent and enzyme treatment	64
12.2.2	Cell lysis by physical disruption.....	65
12.3.	Estimation of total and target protein concentrations.....	65
12.3.1	Express-timator™ Kit protocol.....	65
12.3.2	Estimation of target protein concentration.....	66

Symbol	Description
	Caution, Risk of Danger
	Caution, Risk of Electric Shock
	Protective Conductor Terminal (Ground/Earth)